
## Section 1---Instruction of Assembling Gas Hose

The gas hose can be changed by the customers if there is a need. Please follow the steps below:


1. Mount the Flange Cap to the Gas Hose.

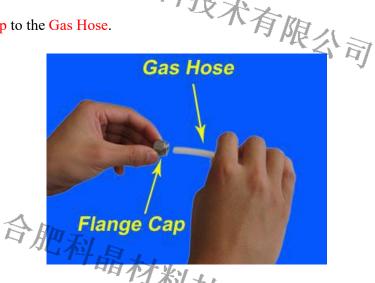


2. Install the Pressing Set onto the gas hose




3. Make the Flange cling to the pressing set.

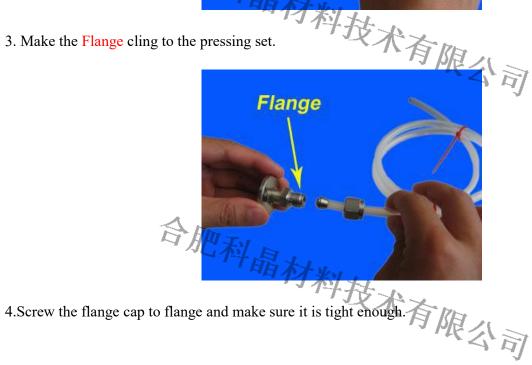

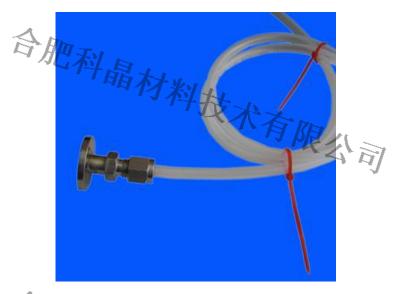

5. Apply epoxy adhesive to glue the Al2O3 ceramic tube and adaptor together.

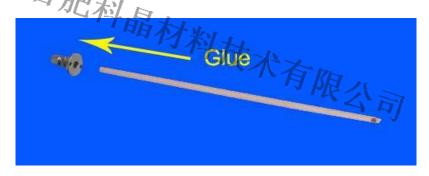


Note: The customer is advised not to change the Al<sub>2</sub>O<sub>3</sub> tube sticking to the flange.

6. Installation is done

Section 2—Instruction of Installing the Gas Purging Adaptor

KJ's muffle furnaces can be easily promoted by installing the gas purging adaptor according to the customers' special use, please follow the steps below to make it right:

1. Unscrew the Two Screws beside the outlet on the top cover of the furnace:

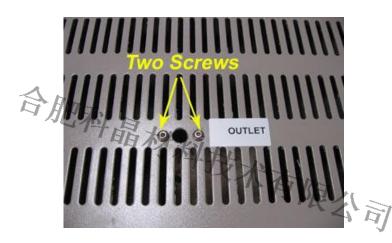


2. Put the Al2O3 Tube labeled "outlet" with the flange and gas hose together into the outlet.



3. Match the screw holes on the flange to the ones beside the outlet, and then tighten the two



4. To install the tube labeled "inlet" to the inlet, repeat the steps 1-3.

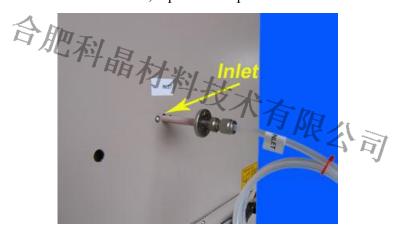

5. Finish the installation and start to use.

